



# AgraQuant® Allergen ELISA Test kits Quick Guide

## AgraQuant® Allergen Test Kits include:

- Set of 5 Standards
- 48/96 antibody-coated microwells
- 1 Bottle of extraction buffer
- 1 Bottle of wash buffer
- 1 Bottle of detection antibody (green cap)
- 1 Bottle of substrate (blue cap)
- 1 Bottle of stop solution (red cap)

## Available kits:

- AgraQuant® Almond 10002021
- AgraQuant® β-Lactoglobulin 10002034
- AgraQuant® Casein 10002037
- AgraQuant® Cashew 10002094
- AgraQuant® Crustacea 10002076
- AgraQuant® Egg white 10002026
- AgraQuant® Fish 10002026
- AgraQuant® Gluten 10002002
- AgraQuant® Gluten G12 10001994
- AgraQuant® Hazelnut 10002006

Shelf life: 12 months

## Test procedures for:

- Raw materials/processed food
- Rinse waters
- Swab samples\*

- AgraQuant® Lupin 10002049
- AgraQuant® Lysozyme 10002090
- AgraQuant® Milk 10002080
- AgraQuant® Mollusk 10006023
- AgraQuant® Mustard 10002070
- AgraQuant® Ovalbumin 10002092
- AgraQuant® Peanut 10001990
- AgraQuant® Pistachio 10002086
- AgraQuant® Sesame 10002064
- AgraQuant® Soy 10002015
- AgraQuant® Walnut 10002030

\* AgraQuant Allergen Swabbing kit (10002205)








## Also Available:

- AgraQuant® Plus Allergen ELISA Test kits
- AgraStrip® Allergens

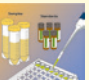



# AgraQuant® Allergen ELISA Test kits Quick Guide

	Almond	β-Lacto-globulin	Cashew	Egg White	Fish	Hazelnut	Lupin	Lysozym	Mustard	Oval-bumin	Peanut	Pistachio	Sesame	Soy	Walnut	Crustacea	Mollusk	Casein	Gluten	Gluten G12	Milk
<b>Extraction</b>	1 g in 20 mL Extraction Buffer at 60 °C															1 g in 20 mL Extraction Buffer (40 °C)	1 g in 20 mL Extraction Buffer (40 °C)	0.5 g in 10 mL Extraction Buffer (60 °C)	1 g in 10 mL 40 % EtOH	0.25 g in 2.5 mL Extraction Solution (50 °C, 40 min) + 7.5 mL 80 % EtOH	0.5 g in 10 mL Extraction Buffer (60 °C)
<b>Sample Dilution</b>	No										1:50		1:10		No (only 100 % meat and sausage samples 1:5)						
<b>Standard Dilution</b>	No										1:100		No		1:100						
<b>1</b>  <b>Add Sample/Standards</b> <b>1<sup>st</sup> Incubation</b>	100 µL Sample/Standard into Well - incubate 20 min in the dark																				
<b>2</b>  <b>Washing</b>	Wash 5x with Wash Buffer																				
<b>3</b>  <b>Add Conjugate</b> <b>2<sup>nd</sup> Incubation</b>	100 µL Conjugate into Well - incubate 20 min in the dark																				
<b>4</b>  <b>Washing</b>	Wash 5x with Wash Buffer																				
<b>5</b>  <b>Add Substrate</b> <b>3<sup>rd</sup> Incubation</b>	100 µL Substrate into Well - incubate 20 min in the dark																				
<b>6</b>  <b>Add Stop Solution</b>	100 µL Stop Solution into Well																				
<b>7</b>  <b>Read with ELISA Reader</b>	Read at 450 nm (with 630 nm differential filter recommended)																				

© Copyright Romer Labs Division Holding GmbH, 2019  
QG\_AQ\_ALL\_EN\_1119\_v01\_MCN